



Widefield 10x eyepiece

Eyepiece tube

Prismatic head

Knurled locking screw, when loosened, permits head to rotate 360° so that eyepiece can be positioned forward or backward for viewing (head retaining screw located on opposite side)

Objective lens 2x

Post locking knob, when loosened, permits microscope to be raised or lowered on post

Focusing knob on both sides of microscope

Support collar stops microscope at lowest point at which microscope can be focused on flat specimen on stage

Stage clips

Stage Plate

Base

Rubber feet prevent sliding, protect table surface